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## BM NA-Off reagent

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Rapid and effective removal of nucleic acid contamination.

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### Usage:

- 1) Clean all kinds of table and instrument surface: Spray the product directly on the table and instrument surface, wipe it with ordinary absorbent paper, and dry it.
- 2) Cleaning of pipette: Disassemble the pipette according to the pipette instruction, soak the part of nucleic acid that is liable to contaminate in the product for 5-10 minutes, then thoroughly soak and rinse with distilled water, dry it, and put it back into the pipette.
- 3) Cleaning of glass and plastic containers: Soak the containers in this product, take them out after 5-10minutes, thoroughly soak and rinse them with distilled water, and then dry them for later use.
- 4) Nucleic acid aerosol cleaning: When there is mild Nucleic acid aerosol pollution in the laboratory, this product can be used to spray the laboratory comprehensively, after 1-2 hours, and with 1:50 diluted 84 disinfectant to wipe the floor and table surface can effectively remove Nucleic acid aerosol pollution; When the laboratory Nucleic acid aerosol pollution is serious, 2-3 times of treatment is recommended, and the treatment effect can be detected after the treatment is completed.